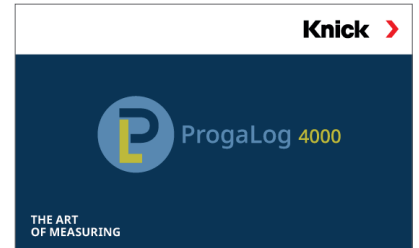


**Headquarters**Beuckestr. 22 • 14163 Berlin  
GermanyPhone: +49 30 80191-0 • Fax: +49 30 80191-200  
info@knick.de • www.knick.de**Local Contacts**

www.knick-international.com

Installationsanleitung  
Installation GuideProgaLog 4000  
PC-Parametriertool für Transmitter  
PC configuration tool for transmitters

Copyright 2021 • Subject to change

This document was last updated on Nov. 18, 2021

The latest documents are available for download on our  
website under the corresponding product description.

099231 TI-201.016-KNXX01

Vor Installation und Verwendung lesen.  
Für künftige Verwendung aufbewahren.  
Read before installation and use.  
Keep for future use.

www.knick.de

**Lizenzschlüssel**

Die PC-Software ProgaLog 4000 startet nach der Installation zunächst im DEMO-Modus.

Mit der Eingabe des Lizenzschlüssels wird der DEMO-Modus beendet. Die  
PC-Software steht vollumfänglich zur Verfügung.**License key**Following installation, the ProgaLog 4000 software initially starts in DEMO  
mode.Entering the license key terminates DEMO mode. The PC software is fully  
available.**Lizenzschlüssel • License key**

## Voraussetzungen

### Hardware

- CPU: 1 GHz Pentium oder vergleichbarer Prozessor
- RAM: 512 MB
- Grafikkarte: 1024 x 768 True Color, 32-Bit
- USB 2.0
- Festplatte: 700 MB

### Betriebssystem:

- Windows 7\*/8/10 (32-Bit- und 64-Bit-Version)
- Microsoft .Net Framework 4.6  
(in Windows 8/10 bereits enthalten)

\*) Bei Verwendung eines PCs mit Windows 7 stellen Sie vor dem Installieren von ProgaLog 4000 sicher, dass Microsoft .Net Framework 4.6 installiert ist (kostenloser Download unter [www.microsoft.com](http://www.microsoft.com))

## Installation von ProgaLog 4000

01. USB-Stick mit einem USB-Anschluss des Computers verbinden.
02. Datei **ProgaLog4000Setup.exe** vom USB-Stick aus ausführen.
03. Installationssprache auswählen.
04. Den Angaben auf dem Bildschirm folgen.
05. Installation fertigstellen.
06. Starten der Software durch Doppelklick auf **ProgaLog-4000Setup.exe**. Beim ersten Mal startet die Software im DEMO-Modus.
07. Zum Beenden des DEMO-Modus Lizenzschlüssel eingeben.

Dokument gültig ab ProgaLog 4000, SW-Version 1.0.0

## System Requirements

### Hardware

CPU: 1 GHz Pentium or comparable processor

- RAM: 512 MB
- Graphics card: 1024 x 768 true color (32-bit)
- USB 2.0
- Hard Disk: 700 MB

### Operating system:

- Windows 7\*/8 or 10 (32-bit or 64-bit version)
- Microsoft .Net Framework 4.6  
(already included in Windows 8/10)

\*) If you have a PC with Windows 7, first make sure that Microsoft .Net Framework 4.6 (free download from [microsoft.com](http://microsoft.com)) is installed before you start the ProgaLog 4000 installation.

## Installing ProgaLog 4000

01. Connect the USB stick to a USB port on the computer.
02. Run the file **ProgaLog4000Setup.exe** from the USB stick.
03. Select the installation language.
04. Follow the instructions on the screen.
05. Complete the installation.
06. Start the software by double-clicking on **ProgaLog4000Setup.exe**. The first time, the software starts in DEMO mode.
07. Enter the license key to terminate DEMO mode.