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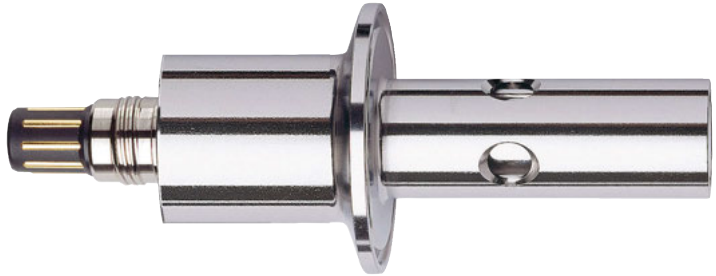
---

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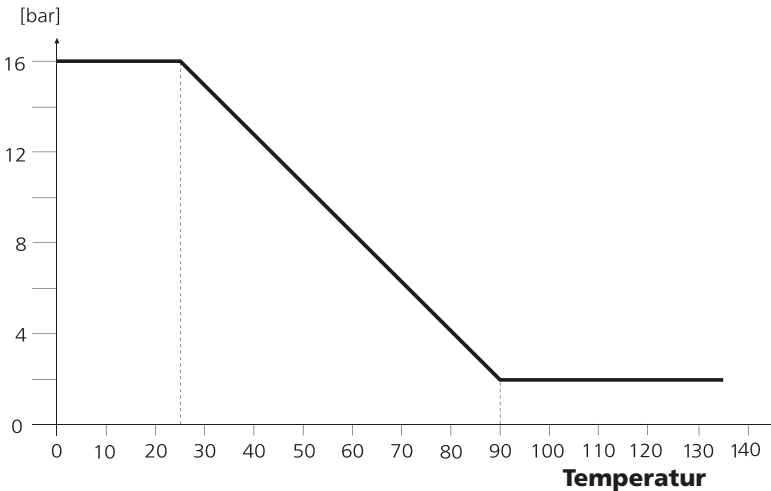
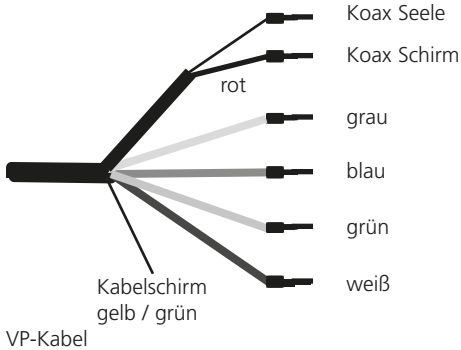
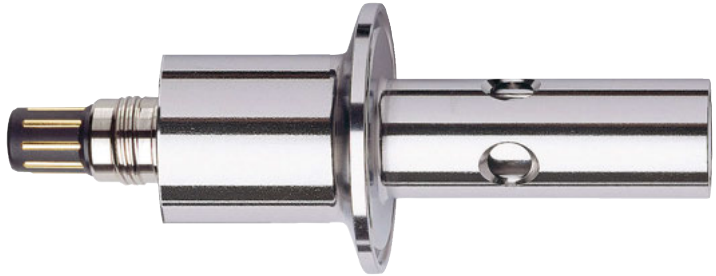
## SE 620 2-Electrode Conductivity Sensor

---

**Be sure to read and observe the following safety instructions!**

### **Warning!**

- When these instructions are not observed, process media under pressure can escape and cause severe injuries.
- Observe correct installation (see dimension drawing)! Otherwise, the assembly is not tight.
- Do not dismantle the sensor as long as the process medium is under pressure!

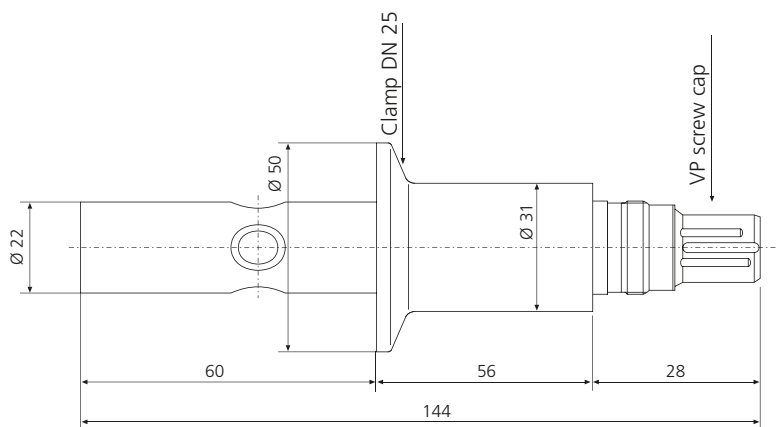
**Knick** ➤

## 1. Applications

The hygienic SE 620 conductivity sensor has coaxially arranged electrodes made of stainless steel. A temperature detector is integrated in the sensor. Electrical connection is made via a VP (VarioPin) cap. Its low surface roughness, the physiologically harmless and FDA-compliant materials, and steam sterilizability make the sensor particularly suitable for the following applications:

Pure and ultrapure water, water for injection (WFI) (pharma), food production, ion exchangers, reverse osmosis, and chip production

## 2. Dimension Drawing

## 3. Calibration / Adjustment

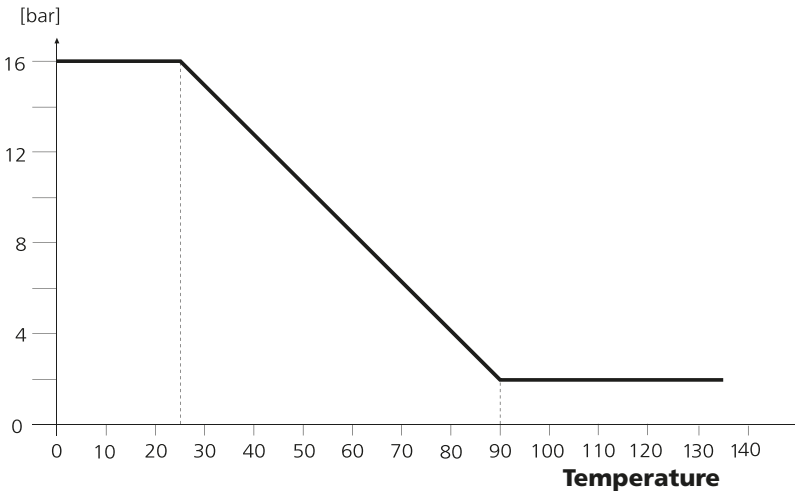
Remove sensor, clean, rinse, and dry it, then immerse it in a calibration solution. Suitable calibration solution ZU 0350 (15  $\mu\text{S}/\text{cm}$ ), ZU 0702 (147  $\mu\text{S}/\text{cm}$ ). Perform a product calibration if the sensor can or shall not be removed. Observe instruction manual of meter. The cell constant of the sensor can be certified by the manufacturer (ZU 0320). For that purpose, you must send in the sensor.

## 4. Specifications

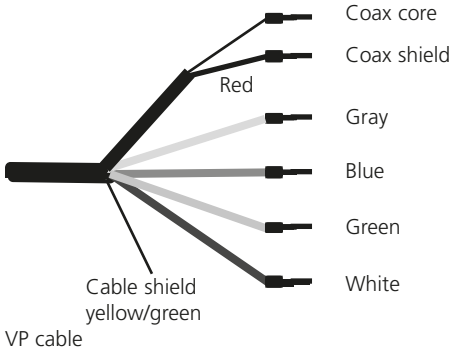
Cell constant:	0.01 /cm	
Range:	0.05 ... 50 $\mu$ S/cm	
Material:	• Sensor incl. electrodes:	Stainless steel 1.4435, electropolished
	• Insulator and O-rings:	FDA-listed plastics
Roughness:	< 0.8 $\mu$ m	
Temperature detector:	Pt 1000	
Temperature:	0 ... 135 °C (steam sterilizable)	
Pressure:	16 bars (25 °C), 9 bars (60 °C)	
Process connection:	Model SE 620/1: Clamp DN 25	
Electrical connection:	VP (VarioPin)	
Dimensions:	See dimension drawing	
Weight:	Approx. 0.5 kg	

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